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1. THE FUNCTION OF KEYS

| | |
|---|---------------------------------|
|  | Turn on/off the balance. |
|  | Change the weighing unit. |
|  | Tare the balance. |
|  | Switch between different modes. |
|  | Enter span calibration. |

2. INSTALLATION

Installs the balance in a firm weighing site and avoid vibration, air flow from air-conditioner, open window, or ventilator, and electromagnetic waves.

Note: Before installation and use: please make sure to loosen the red screw which is use to protect the transducer. And please keep the red screw properly in order to use it when the balance need to transport.

3. TURNING ON THE POWER

Insert the plug of the AC adapter into the DC IN connector on the rear of the balance or dry batteries of 12v into the battery compartment.

Turn on the  key, and it will display “XXXX” which stands for the maximum weighing value, and wait for seconds (depending on the working environment) to display “0--0.0--0.00--0.000” weighing mode.

4. SPAN CALIBRATION

4.1 NOTA

Please allow the balance to warm up for 10-30 minutes (depending on the different type of balances) when first time turn on  in order to make the balance adapt the environment.

4.2. CALIBRATION

- a. Verify that the pan is empty, press and hold  key about 5 seconds, when the display shows “CAL”, release the  key.
- b. Load the indicated weight which's number is blinking on the LCD, wait the number stop blinking, Un-

load the weight.

- c. "---" is displayed briefly to indicate completion of span calibration and soon the balance returns to waiting mode.
- d. Repeat calibration several times, Do it as slowly as possible or operate linear calibration after first calibration.

4.3. LINEAR CALIBRATION

- a. Verify that the pan is empty, press and hold  key about 5 seconds, when the display shows "CAL", release the  key.
- b. Press and hold  key about 5 seconds, the display starts to show the every session of linear calibration values, the blinking number indicates which weight should be placed on the pan.
- c. Load the indicated weight, and wait for the display number stops blinking, it means the balance has record this linear value. Unload the weight.
- d. Wait for the next blinking number, repeat step "c" to calibrate the next value.
- e. When the display briefly shows "0.000", it means the linear calibration is completed.

5. BALANCE COUNTING OPERATION

5.1 BALANCE COUNTING

- a. Verify the balance is on the waiting mode (tare) which the display shows "0.000", and shortly press  key, the display will show "10" and blink. The "g" on the right-bottom will change to "pcs".
- b. Press the  key to select a sample size which will change sequentially as following 10-20-50-100-150-200-250-500.
- c. Place the same number of samples you selected on the pan then press  key into the counting operation.

5.2 NOTA

User can press  key again to exit the counting function and return to the waiting mode.

6. CHANGE THE UNIT DISPLAY

Press the  key to changes it as following sequence "CT" (carat), "OZ" (ounce), "1b" (pound), "g" (gram). The default unit is "g".

7. SPECIFICATION

| MODEL | ABS6 | ABS30 |
|-------------------|-----------------|-------|
| Capacity (kg) | 6 | 30 |
| Fraction (g) | 0,05 | 0,1 |
| Pan size (mm) | 2115 x 295 | |
| Dimensions (mm) | 310 x 310 x 115 | |
| Net Weight (kg) | 3,9/5 | |
| Warm up (minutes) | 10-20 | |

8. SAFETY PRECAUTIONS

Allow the balance to Warm up and calibrate before starting to use the balance. Use the calibration weight to calibrate the balance with a weighing error. Do not use the balance in hazardous areas. This includes areas where the balance is exposed to dust or flammable gases and liquids.

Troubleshooting

| SYMPTOM | PROBABLE CAUSE(S) | COUNTERMEASURE |
|------------------------------------|--|---|
| "---" is displayed on top | The load is too heavy | Use balance within its capacity or re-calibration |
| Display fluctuates | Affected by vibration or air flow | Install the balance in a proper site |
| | You don't loosen the red screw | Loosen the red screw |
| | Windbreak glass window isn't closed | Close the window |
| | The Pan touches the shell or some unexpected item is on the pan. | Check the pan and clean it |
| The weighed result is not accurate | The weighing item is fluctuates | Fix the item |
| | Span calibration has not been done | Calibrate it properly |
| | Tare has not been done | Tare before weighing |
| | The balance is not in the level | Level the balance |

9. STANDARD ACCESSORIES

In addition to a balance, this package includes the following items:

| | |
|-----------------------|--------|
| AC adapter | 1 |
| Pan | 1 |
| Weight | 0- 1-2 |
| Printed documentation | 1 |
| Warranty document | 1 |
| Qualification | 1 |

10. WARRANTY

This scale is warranted against defects of manufacturing and materials for a period of 1 year, from the delivery date.

During this period, GIROPÈS SL will take charge of repairing the scale.

This warranty does not cover defects or damaged caused by misuse, overloading or improper installation contrary to the recommendations described in this manual.

This warranty does not cover shipping costs for the reparation of the balance.

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Declare under our responsibility that the denominated product:

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To wich declaración referes to, conform wich the followings standards or other normative documents:

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2004/108CE directive on the electromagnetic compatibility.

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And the following harmonized norms/ normes harmonisées:

EN61000-6-3: 2007

EN61000-6-1: 2007

EN61000-3-2: 2006+A2: 2009

EN61000-3-3: 2008

Jordi ribalta

Director General
General Manager

Baxtran

marca propiedad de/
est une marque de/
trade mark propriety of:

Pol. Empordà Internacional
Calle F. Parcela 15-16
17469 VILAMALLA
(Girona) SPAIN
T. (34) 972 527 212
F. (34) 972 527 211